# Mini-Lab

## Instruction Manual

2025

Mini-Lab

dupal<sup>UK</sup>

## Instruction Manual

<b>Mini Lab's content</b>	<b>P.3</b>
<b>Incubator instructions</b>	<b>P.4 to 6</b>
- Operation mode	p.4 to 5
- Alarm patterns & recommended actions	p.6

## Notice

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# C4Hydro Mini-Lab Instruction Manual

## The C4Hydro Mini-Lab contains:

Incubator +  
power adapter with plug

50 mL syringe

10 mL syringe with a connector

1 mL syringe with a connector

Cap

### Note:

In order to test water, the Mini-Lab **has to be used with a C4Hydro Water Test**. The content of the Mini-Lab **will be re-used for every test** and should not be discarded after water testing.

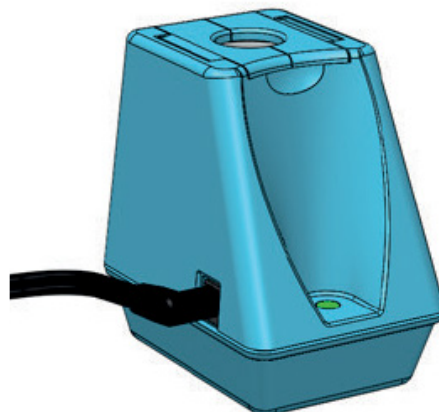
# Incubator instruction manual

The purpose of the incubator is to maintain the contents of a syringe at a constant temperature.

## Operation mode:

- 1) Connect the power adapter to the connector of the incubator. Make sure the 2 half lids are closed.


- 2) Connect the power adapter to a main power supply (AC 100–240V / 50–60Hz).

- The LED light blinks once red, green and blue to confirm power supply.
- The LED light blinks green until the incubation temperature is reached
- The LED light remains green when the right incubation temperature is reached.

- 3) When the LED remains green, open the 2 half lids at the top of the incubator.

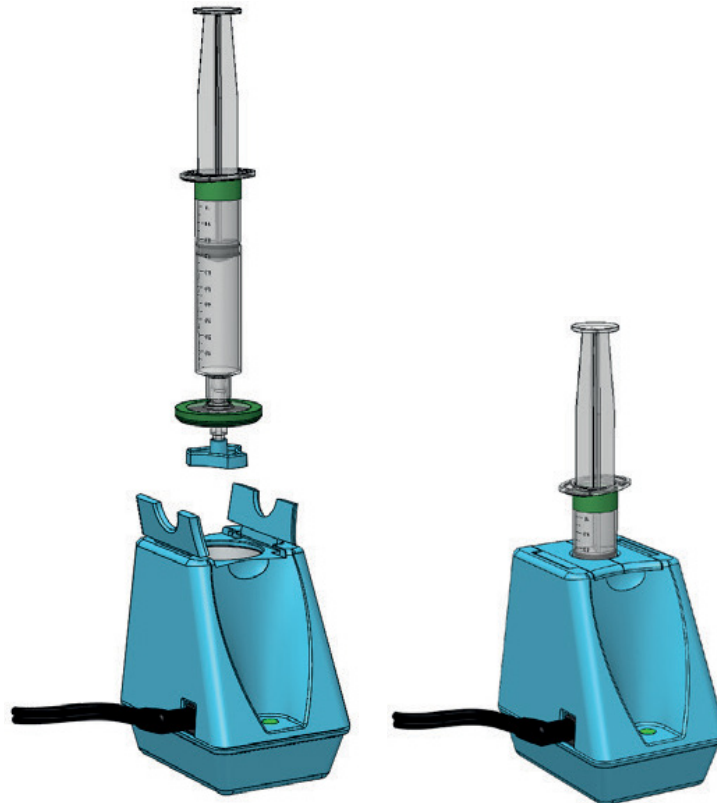- 4) Place the syringe with filter and cap inside the incubator in vertical position as shown.

- 5) Close the 2 half lids at the top of the incubator and do not disconnect the incubator from power supply during incubation.
- 6) When the incubation step is over, open the 2 half lids at the top of the incubator before taking the syringe out.
- 7) When the water sample analysis is over, disconnect the power adapter from the power supply and then disconnect the incubator from the power adapter.

# Incubator instruction manual

The following table describes the alarm patterns and recommended actions to be taken:

INCUBATOR LED COLOR	ALARM MEANING	ACTION
LED LIGHT BLINKS BLUE 	The incubation temperature is not reached after 30 minutes	Ambient temperature of the room is too cold. The incubator should be used in a warmer ambient temperature room.
LED LIGHT BLINKS RED 	Functional issue. The incubator will not heat.	The incubator and/or the adapter should be replaced.

**Electrical specifications:**  
Plug power adapter  
AC 100–240V to DC 12V–2A

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Dupal UK Limited